

EN

EP.V.36622.00 | 05/2025

asecos®



ION_{LINE}

CORE

USER MANUAL

Safety storage cabinets for storage and charging of lithium ion batteries

ION-CORE-90

IO90.195.060.CC.WDC



IO90.195.120.CS.WDC



IO90.195.060.CS.WDC

EN: Content

1. NOTES • GUIDELINES • GUARANTEE	4
1.1. General safety notes	4
1.2. Guarantee	4
1.3. Cabinet Details	4
2. TRANSPORT	5
2.1. Tilting the cabinet	5
2.2. Dismantling of the transport packaging	5
2.3. In-plant transport	5
2.4. Tilting onto the side wall	6
3. INSTALLATION	6
3.1. Alignment Of The Cabinets	6
4. Commissioning	6
4.1. Connection to the power supply	6
4.2. Potential-free alarm contact	7
5. CLOSING	7
5.1. In general	7
5.2. Locker System	8
6. INTERIOR FITTINGS	8
6.1. Bottom collecting sump	8
6.2. Insolated Shelves (height-adjustable)	8
6.3. Total power rating of the power socket strips	9
7. STORAGE	9
7.1. General information on batteries	9
7.2. Notes on storage and charging	9
8. SMOKE DETECTOR • PRESSURE RELIEF	9
8.1. Smoke detector	9
8.2. Pressure relief	10
9. ALARM OVERVIEW	10
9.1. Error and alarm overview	10
10. BATTERY FIRE • EVENT OF FIRE • DISPOSAL	10
10.1. Fire inside the cabinet (battery fire)	10
10.2. Opening the cabinet after the fire	11
10.3. Disposal	11
11. SAFETY CHECKS	11
11.1. Smoke detector maintenance	11
11.2. Cleaning	12
11.3. Contact	12
12. TECHNICAL DATA	12
13. TECHNICAL DRAWING	12

DE: Inhalt

1. HINWEISE • RICHTLINIEN • GEWÄHRLEISTUNG	13
1.1. Allgemeine sicherheitstechnische Hinweise	13
1.2. Gewährleistung	13
1.3. Schrankdetails	13
2. TRANSPORT	14
2.1. Umlegen des Schrankes	14
2.2. Demontage Transportverpackung	14
2.3. Innerbetrieblicher Transport	14
2.4. Auf Seitenwand kippen	15
3. AUFSTELLUNG	15
3.1. Ausrichten der Schränke	15
4. INBETRIEBNAHME	15
4.1. Anschluss an die Stromversorgung	16
4.2. Potentialfreier Schaltkontakt	16
5. SCHLIESSUNG	17
5.1. Allgemein	17
5.2. Schliessfachsystem (optional)	17
6. INNENAUSSTATTUNG	17
6.1. Bodenauffangwanne	17
6.2. Isolierte Fachböden (höhenverstellbar)	18
6.3. Gesamtleistung der Steckdosenleisten	18
7. LAGERUNG	19
7.1. Allgemeine Hinweise zu Batterien	19
7.2. Hinweise zu Lagerung und Laden	19
8. RAUCHMELDER • DRUCKENTLASTUNG	19
8.1. Rauchmelder	19
8.2. Druckentlastung	19
9. ALARMÜBERSICHT	20
9.1. Störung und Alarmübersicht	20
10. AKKUBRAND • BRANDFALL • ENTSORGUNG	20
10.1. Brand im Schrank	20
10.2. Öffnen des Schrankes nach dem Brand	20
10.3. Entsorgung	21
11. SICHERHEITSTECHNISCHE ÜBERPRÜFUNG	21
11.1. Rauchmelderprüfung	21
11.2. Reinigung	21
11.3. Kontakt	22
12. TECHNISCHE DATEN	22
13. TECHNISCHE ZEICHNUNG	22



asecos GmbH
Customer service
Weierfeldsiedlung 16–18
D-63584 Gründau

Fax: +49 60 51 - 92 20-10
email: service asecos.com

YOUR PERSONAL DOCUMENTATION TO THE asecos SAFETY CABINET

Dear Customer,

you have made a decisive investment in safety for your company by purchasing this asecos safety storage cabinet. You now own an innovative product made of top-quality materials guaranteeing the highest quality standards.

asecos safety storage cabinets have complete authorisation documents. We archive the authorisation documents for every individual cabinet, keeping them ready for you should you ever need them (e.g. for a works inspection or similar). Simply request them using this form.

Tear of or copy that page and return to us by fax with your address and serial number of the cabinet on it.

Yours sincerely
asecos GmbH

Contact

Company		
<input type="text"/>		
Street	Postal code	Town
<input type="text"/>	<input type="text"/>	<input type="text"/>
Name of contact person		
<input type="text"/>		
email	Phone No.	
<input type="text"/>	<input type="text"/>	
Serial numbers of safety storage cabinets		
<input type="text"/>	<input type="text"/>	<input type="text"/>
<input type="text"/>	<input type="text"/>	<input type="text"/>
<input type="text"/>	<input type="text"/>	<input type="text"/>
<input type="text"/>	<input type="text"/>	<input type="text"/>

OPERATING INSTRUCTIONS

Dear customer,

Thank you very much for purchasing a safety storage cabinet from our company, with which you have made a decisive investment in the safety within your company. Our safety storage cabinets make the storage of hazardous materials at the workplace safe and convenient for you.

Please read these operating instructions very carefully. Get to know the advantages and simple operability of our safety storage cabinets in detail. This simplifies the daily handling of hazardous materials for you.

Many thanks
Your asecos team


- 1. NOTES • GUIDELINES • GUARANTEE5**
 - 1.1. General safety notes 5
 - 1.2. Safety instructions for the storage of lithium-ion batteries 5
 - 1.3. Guarantee 5
 - 1.4. Cabinet Details 5
- 2. TRANSPORT6**
 - 2.1. Tilting the cabinet 6
 - 2.2. Dismantling of the transport packaging 6
 - 2.3. In-plant transport 7
 - 2.4. Titling onto the side wall 7
- 3. INSTALLATION7**
 - 3.1. Alignment Of The Cabinets 7
- 4. COMMISSIONING7**
 - 4.1. Connection to the power supply 8
 - 4.2. Potential-free alarm contact 8
- 5. CLOSING9**
 - 5.1. In general 9
 - 5.2. Locker System 9
- 6. INTERIOR FITTINGS9**
 - 6.1. Bottom collecting sump 9
 - 6.2. Insolated Shelves (height-adjustable) 9
 - 6.3. Total power rating of the power socket strips 10
- 7. STORAGE10**
 - 7.1. General information on batteries 10
 - 7.2. Notes on storage and charging 10
- 8. SMOKE DETECTOR • PRESSURE RELIEF11**
 - 8.1. Smoke detector11
 - 8.2. Pressure relief11
- 9. ALARM OVERVIEW11**
 - 9.1. Error and alarm overview11
- 10. BATTERY FIRE • EVENT OF FIRE • DISPOSAL11**
 - 10.1. Fire inside the cabinet (battery fire)11
 - 10.2. Opening the cabinet after the fire 12
 - 10.3. Disposal 12
- 11. SAFETY CHECKS12**
 - 11.1. Smoke detector maintenance 12
 - 11.2. Cleaning 13
 - 11.3. Contact 13
- 12. TECHNICAL DATA13**
- 13. TECHNICAL DRAWING14**

1. NOTES • GUIDELINES • GUARANTEE

1.1. GENERAL SAFETY NOTES

- When handling lithium-ion batteries, observe the applicable regulations and the information in these operating instructions
- Work on the electrical system is to be carried out only with the power turned off and only by qualified electricians – refer here to the regulations of the local electricity supply company.
- General damage to electronic components is to be repaired without delay by an asecos employee.
- Use only intact and undamaged mains cables for the battery charger
- Electrical protection in accordance with local standards must be provided by the customer (cabinets do not have their own RCD circuit breaker or circuit breaker)
- The on-site installation conditions are to be observed.
- The instructions of the supervisory engineering department must be followed.
- Observe accident prevention regulations and workplace ordinance
- **Ensure that the necessary safety checks are only carried out by authorised staff using original spare parts**
- Only use the cabinet after having been properly instructed; access is to be forbidden to unauthorised persons.
- The doors are permanently self-closing and must not be pushed shut manually
- The pivoting area of the doors is to be kept free at all times; doors are to be kept closed
- By assigning trained/authorised technical personnel you can prevent the malfunctions, damage and corrosion damage that result from inappropriate transport.
- Observe the upper limits for stored quantities, loading etc.
- The following substances may not be stored in the cabinets with a fire suppression system: acids, alkalis, magnesium, other metals (in powder form)
- If cable openings are no longer required due to new usage circumstances, they must be closed in accordance with the manufacturer's specifications.
- **Observe the notes on maximum size and general storage of the batteries in these instructions**

Set-up requirements

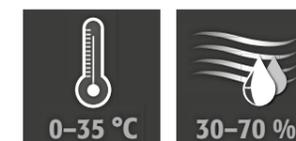

1.2. SAFETY INSTRUCTIONS FOR THE STORAGE OF LITHIUM-ION BATTERIES

- Any charger included will switch off in the event of fire or overheating
- Follow the instructions for storing your batteries in accordance with the battery manufacturer's operating instructions
- Batteries can produce toxic gases, therefore connection to an exhaust air system is recommended.
- Alternatively, a risk analysis must be carried out. The result determines whether additional safety measures are necessary to minimize the risk of toxic, flammable gases in the work area.

1.3. GUARANTEE

The guarantee for this product is agreed between you (the customer) and your dealer (the seller). As the manufacturer, asecos guarantees the products listed in the operating instructions for a period of 24 months from the date of delivery. All model safety equipment are subject to a compulsory annual inspection by specialised staff authorised by the manufacturer. Otherwise the customer's guarantee claim against the manufacturer expires. Please note that the warranty is also void if holes are drilled or modifications are made without consulting the manufacturer asecos.

1.4. CABINET DETAILS

Cabinet data: logbook (included with the cabinet)
Technical drawing: see appendix
Technical data: table in appendix

ION-CORE-90

Safety storage cabinets for lithium-ion batteries
Comprehensive fire protection with the proven evacuation and alarm forwarding concept. During active storage, lithium-ion batteries or battery packs are charged or partially discharged (60-70%) in the cabinet using a charger.

Model	Lithium ion batteries		Integrated technical ventilation	Fire suppression unit	Alarm system
	Storing	Charging			
IO90.195.060.CC.WDC	✓	✓			✓
IO90.195.120.CS.WDC	✓				✓
IO90.195.060.CS.WDC	✓				✓

2. TRANSPORT



CAUTION:

Transport the cabinet in an upright position on a pallet truck, tied and secured against slipping, until the final place of installation is reached. The transport locks in the door joints may only be removed directly at the place of installation! Inappropriate transport can lead to concealed damage to the fire protection insulation! We can only guarantee the necessary quality if the cabinet is transported to the place of its use by our specially trained staff.


2.1. TILTING THE CABINET



CAUTION:

Tilting the cabinet may only be done without jolts! When using a stair robot, observe the maximum load capacity of the device according to the manufacturer's instructions!

2.2. DISMANTLING OF THE TRANSPORT PACKAGING

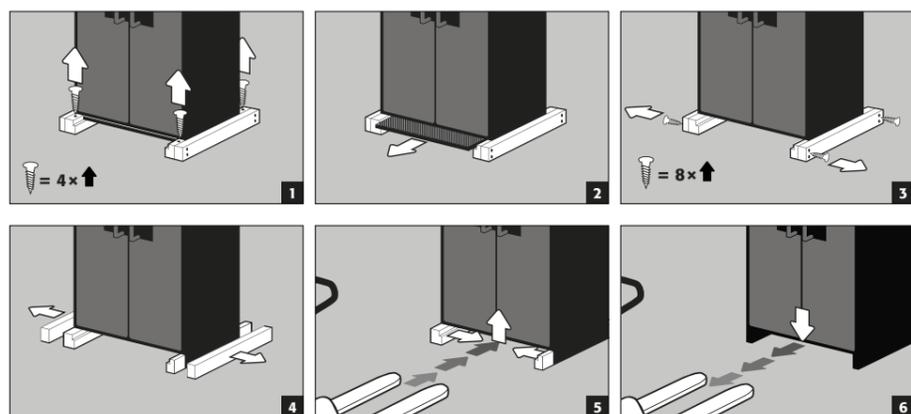


CAUTION:

Cabinets with a width of 600 mm: The clear entry width of the base is 520 mm. Please note this when choosing your pallet truck! Devices with widths greater than the entry widths must not be used.


2.3. IN-PLANT TRANSPORT

- In-plant transport is also possible without transport locks (inserted as standard in the door joints)

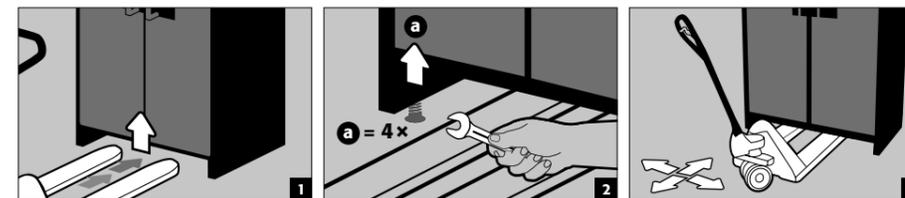


2.4. TITLING ONTO THE SIDE WALL

- Tilting onto the side wall is only possible with the optionally available tilting bracket (order no. 29556).

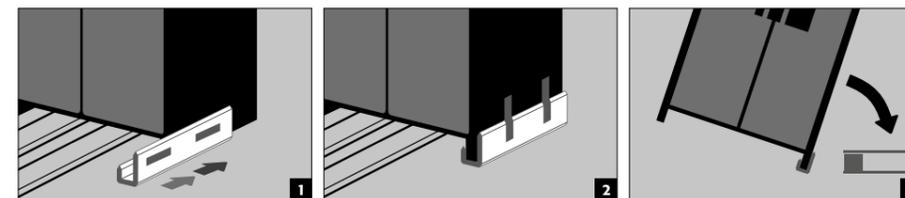


3. INSTALLATION

3.1. ALIGNMENT OF THE CABINETS

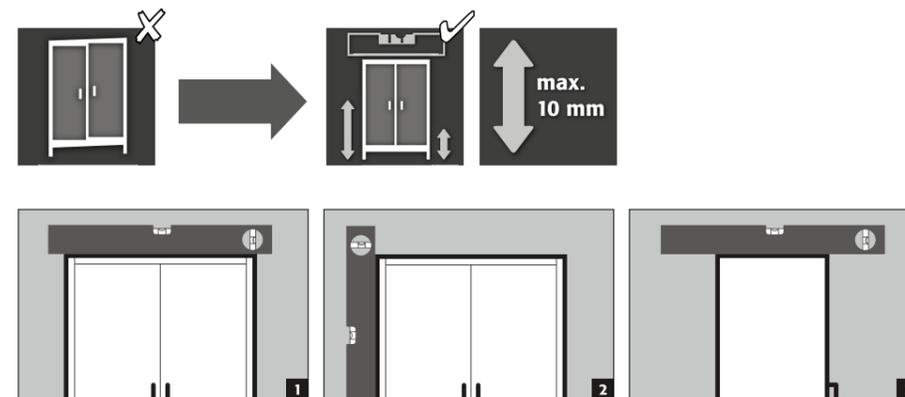

CAUTION:


Door elements must not scrape against the fire prevention seals in the fold of the door when opening and closing! Doors with an automatic closing mechanism must close automatically from every position and the lock must be able to lock!

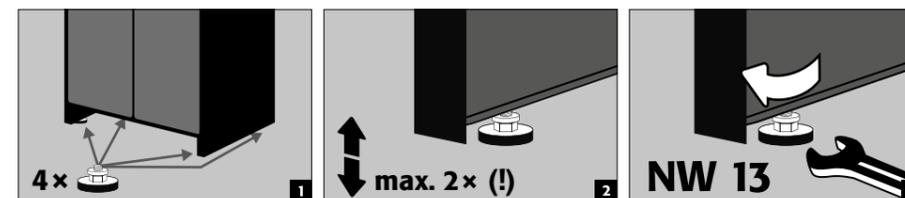


4. COMMISSIONING

- Before putting into operation for the first time, the user must carry out an examination of the safety storage cabinet for possible damage, such as defective or loose sealing elements, correct alignment and perfect functioning of the door elements. Use the cabinet and accessories only if they are in an orderly condition.


4.1. CONNECTION TO THE POWER SUPPLY

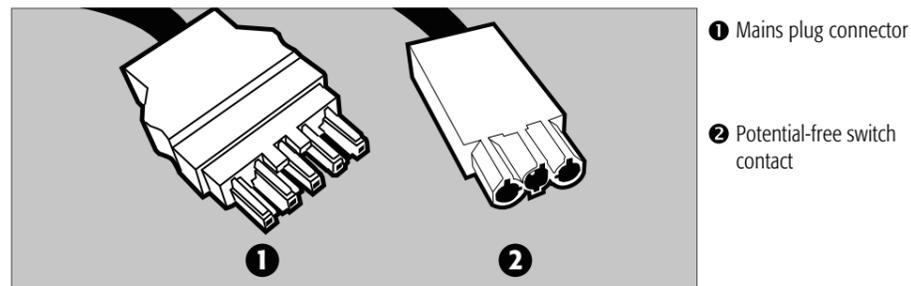

Connection to the power supply

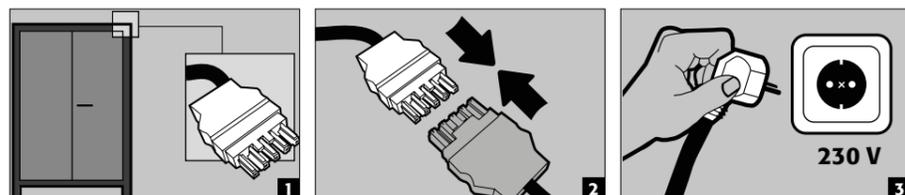

Connection to the power supply with 400 V (optional with item 38038)

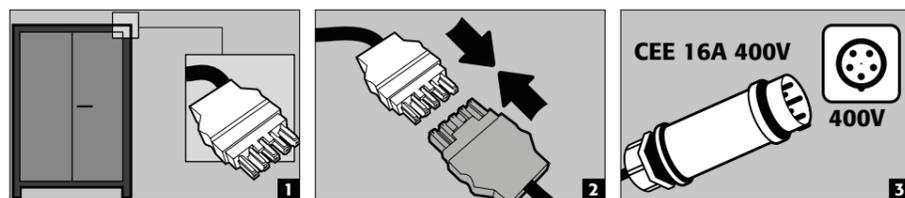


NOTE:
Retrofitting is easy due to the plug-in connection, so that no intervention in the electrical components is necessary. The power supply must be fused on site with a maximum of 16A. Electrical protection in accordance with local standards must be provided on site. The cabinets do not have their own RCD circuit breaker or circuit breaker.

4.2. POTENTIAL-FREE ALARM CONTACT



NOTE:
The potential-free alarm contact is used to connect a signal to a control centre/control room. Direct integration into a fire alarm control panel (FACP) is not recommended or may only be carried out in consultation with the person responsible for the system.
However, it is always recommended to connect the signal to a manned control centre/control room!
The potential-free switching contact must always be connected by the customer (not a service).

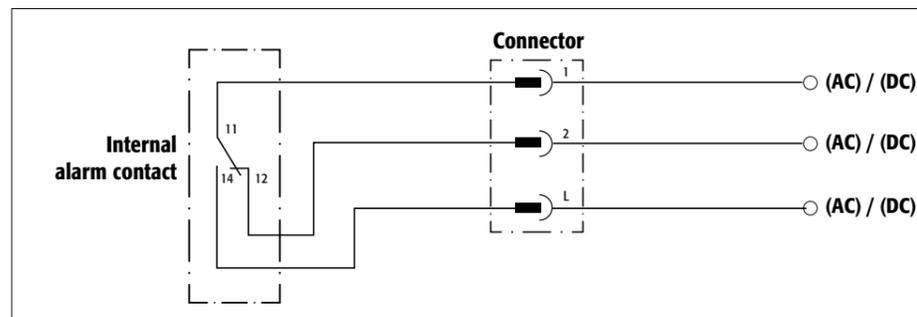


Connection instructions

- Use only the supplied mating part (colour-coded brown) to the plug for the connection.
- The connection must be done by a qualified electrician.
- The internal switch contact is designed for a max. DC voltage of 24 V or a max. AC voltage of 230 V.
- The maximum current load is 5 A at 230 V AC and 10 A at 24 V DC.
- The internal switch contact is a changeover contact; in case of alarm, therefore, the switching state may be queried as „opened“ or „closed“.

5. CLOSING

5.1. IN GENERAL

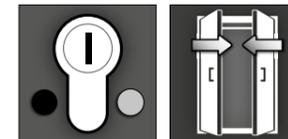

The doors are permanently self-closing and self-locking. They must not be closed manually. The cabinets have a profile cylinder lock with locking status indicator and can be integrated into a locking system. A profile half cylinder (30/10) with adjustable cam must be used.



ATTENTION:
The owner/user must ensure that all doors are kept closed whenever the contents of the cabinet are not being accessed. In general, it must be noted that the cabinets do not possess an emergency unlocking facility. This means that persons trapped inside the cabinet cannot free themselves!

5.2. LOCKER SYSTEM

- Lockers can be closed manually and each has a cylinder lock with its own key pair
- An additional master key opens all 7 lockers
- Lockers and keys can be individually numbered using the key ring set and sticker sheet provided

6. INTERIOR FITTINGS


6.1. BOTTOM COLLECTING SUMP

Leaks:

- Liquid in the sump is to be collected using suitable means.
- The choice of means is your own responsibility.



6.2. INSOLATED SHELVES (HEIGHT-ADJUSTABLE)

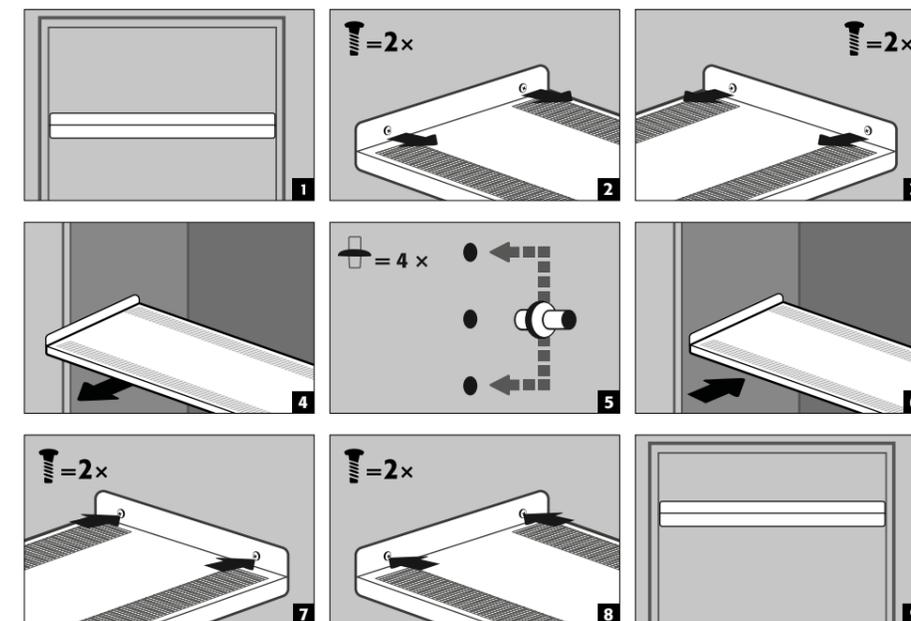

LOAD CAPACITY (KG/ DRUM PLATFORM)

IO90.195.120.CS.WDC	IO90.195.060.CC.WDC IO90.195.060.CS.WDC
 max. 75 kg	 max. 25 kg



CAUTION:
Please note that dynamic forces act when loading the cabinets. Always place batteries carefully in the cabinet!



CAUTION:
The position of the shelves / 2nd level drawer and socket strips cannot be changed.

6.3. TOTAL POWER RATING OF THE POWER SOCKET STRIPS

Standard: single-phase, 230 V

Version	EU	CH	UK	FR	other regions:
fusing	16 A	10 A	13 A	16 A	Please get in touch with your asecos contact person. The maximum power and protection may differ here.
max. total power	3,68 kW	2,3 kW	2,99 kW	3,68 kW	

Optional: 2-phase, 400 V (accessories article 38038)

Version	EU	CH	UK	FR	other regions:
fusing	2 x 16 A	2 x 10 A	2 x 13 A	2 x 16 A	Please get in touch with your asecos contact person. The maximum power and protection may differ here.
max. total power	7,36 kW	4,6 kW	5,98 kW	7,36 kW	



ATTENTION:
The load on the system is to be distributed as evenly as possible over the power socket strips! The individual power strip must not be loaded with more than the specified power max. (see table)!
The necessary fuse protection is to be provided by the customer!

7. STORAGE

7.1. GENERAL INFORMATION ON BATTERIES



ATTENTION:
Never store obviously damaged lithium-ion batteries inside buildings. Dispose of them without delay in disposal containers that are provided outside the building and approved for transport.



CAUTION
Only batteries with a maximum weight of 15 kg may be stored in the cabinets.

7.2. NOTES ON STORAGE AND CHARGING

Storage

- It is recommended to store new and used lithium-ion batteries separately (each on a different storage level) in the safety storage cabinet.

Occupation of the storage levels

For this basic version, asecos recommends a load of 1 kWh per storage level.

8. SMOKE DETECTOR - PRESSURE RELIEF

8.1. SMOKE DETECTOR

These models have a battery-operated smoke detector.

Battery change



CAUTION:
The use of rechargeable batteries is not permitted!
The life of the battery is highly dependent on local conditions such as temperature, temperature fluctuations, humidity and the number of test alarms/alarms, among other things. In the case of lithium, this is up to 5 years. The smoke detector announces a necessary battery change approx. 30 days in advance (see 10.2).



8.2. PRESSURE RELIEF

All ION-LINE models have a pressure relief flap built into the head section, which opens briefly in the event of a pressure increase and thus releases the pressure from the cabinet.

9. ALARM OVERVIEW

9.1. ERROR AND ALARM OVERVIEW

EVENT	LED RED ON SMOKE DETECTOR	ACOUSTIC ALARM	ACTIONS
Smoke detector detects smoke in the cabinet	flashing	pulsating alarm tone	see 12.1
Triggered by connected detectors	off	pulsating alarm tone	The triggering detector can be identified by parallel to the alarm tone flashing LED
Battery replacement due	flashing	short beep every 45 seconds	see 8.3
Operational readiness	flashes every 45 seconds	off	
Malfunction	flashes alternately with the beep	short beep every 45 seconds	change smoke detector

10. BATTERY FIRE - EVENT OF FIRE - DISPOSAL

10.1. FIRE INSIDE THE CABINET (BATTERY FIRE)



- For fast transport the cabinets are equipped with a transport base. The cabinets are automatically disconnected from the mains supply in the case of transport.




NOTE for tall cabinets

Evacuation by at least 2 persons is recommended. Only qualified personnel (e.g. fire department) may carry out transportation in the event of a fire.



CAUTION:
The doors must be locked before transport! Depending on the door heights, it may be necessary to remove the ventilation attachment beforehand. Transport may only be carried out by qualified personnel!

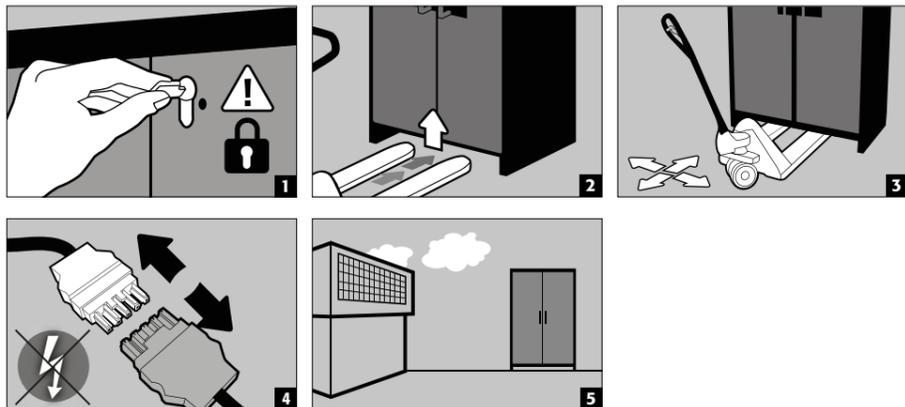

10.2. OPENING THE CABINET AFTER THE FIRE



CAUTION:
Do not open the cabinet until it has cooled down. This is 6 times the fire duration! The cabinet may only be opened by authorised personnel (e.g. fire brigade)!
Depending on the duration of the fire, an ignitable vapour-air mixture may have formed, therefore remove all ignition sources within a 10-metre radius around the cabinets before opening. Use only non-sparking tools! Open the cabinets with extreme caution!

10.3. DISPOSAL



The models can be disposed of once they have been dismantled and the materials sorted.


11. SAFETY CHECKS

As safety equipment the cabinets have to be checked for safety at least once per year. The next checking date can be taken from the service sticker on the outside of the door. This annual check can be carried out with the necessary care, and for securing your warranty claims in the case of fire, only by an authorised asecos employee (refer also to our service brochure regarding this).
A necessary service is indicated by service sticker on the door of the cabinet. Within the context of the annual check, all safety-related parts, the smoke detector and the alarm signalling will be checked.

11.1. SMOKE DETECTOR MAINTENANCE



ATTENTION
In accordance with DIN 14676 the proper function of the smoke detector must be checked at least once per year.

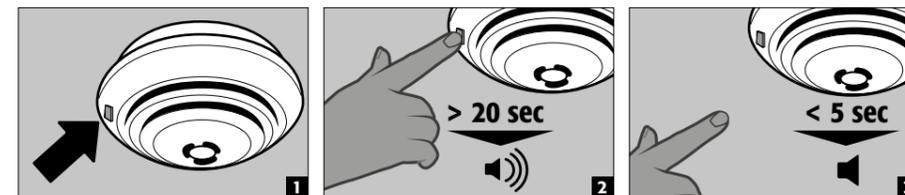

- The smoke detector is completely tested with the LED test button (fig. 1): battery function test, electronic smoke chamber test and a test of the evaluation electronics.
- After releasing, the test alarm resets itself
- Following a successful test, the alarm is silenced and the LED flashes every 45 seconds - the smoke detector is ready to operate
- If the test failed, see error and alarm overview for error analysis

Self-test

- The smoke detector carries out an automatic self-test, in which the evaluation electronics as well as the voltage and the internal resistance of the battery are checked about every 45 seconds.
- This check is signalled by a short flashing signal of the red LED.

11.2. CLEANING

The cabinets can be cleaned with a mild household cleaner and a soft cloth. In case of damage please contact your dealer in order to have the cabinet repaired using original spare parts.

11.3. CONTACT



CONTACT:
In the case of defects or complaints about our products (within and also after the warranty period), and for requesting safety checks or taking out a service contract, please contact our service hotline on:
Tel: +44 1785 22 70-90 | info@asecos.co.uk (for great Britain and Ireland)
Tel: +49 1805 92 20 92 | service@asecos.com (international)

12. TECHNICAL DATA

ION-CORE-90		IO90.195.060.CC.WDC	IO90.195.060.CS.WDC	IO90.195.120.CS.WDC
Type		90	90	90
External dimensions W x D x H	mm	599 x 615 x 1953	599 x 615 x 1953	1193 x 615 x 1953
Internal dimensions W x D x H	mm	450 x 503 x 1647	450 x 503 x 1647	1050 x 503 x 1647
Weight without interior equipment	kg	305	300	480
Distributed load	kg/m ²	935.50	930.00	560.55
extraction air connection	DN	75	75	75
Entry width transport base	mm	526	526	1120
Entry height transport base	mm	90	90	90
Max. shelf load (evenly distributed)	kg	25	25	75


Electrotechnical data

Nominal voltage	V	230/400
Frequency	Hz	50/60


Total power rating of the power socket strips

		EU	CH	UK	FR/BE
Fuse (1-phase)	A	16	10	13	16
Power max. (1-phase)	kW	3,68	2,3	2,99	3,68
Fuse (3-phase)	A	2 x 16	2 x 10	2 x 13	2 x 16
Power max. (3-phase)	kW	7,36	4,6	5,98	7,36


13. TECHNICAL DRAWING

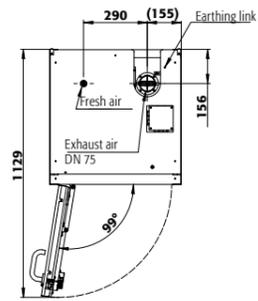

IO90.195.060.CC.WDC

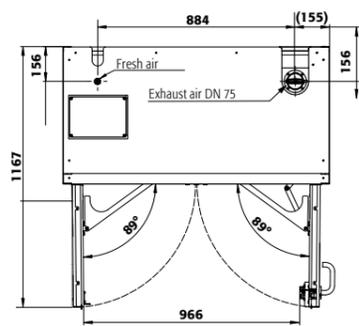

IO90.195.120.CS.WDC

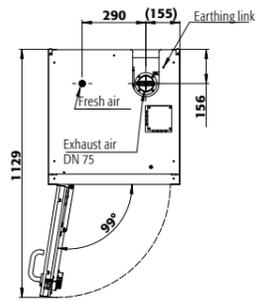

IO90.195.060.CS.WDC

